

NIFTILIDS™

- Top quality specimen containers made from opaque polypropylene
- British made
- Leak-proof lids to meet COSHH regulations
- Lids easy to use and usable time after time
- Colour coded lids available

Please see inside for further details.



HOW TO



OPENING

When removing the lids, it is important that the lip is folded back **all the way round** for the lid to then effortlessly lift off. Failure to do this by tugging, tearing or pulling the lid off will damage the lid and rim of the container, often weakening the flange and losing its water-tight seal.



CLOSING WITH LIP TURNED DOWN (for the first time)

When sealing the containers with the lid, ensure that the rim is pushed down firmly. Then to complete the watertight seal, place the container on a level that you are able to lean over it and using your upper body weight through your arms place your hands side by side with thumbs touching and press down hard on the rim in a sharp movement with the balls of your hands until a “snap” is heard. This forms the watertight seal. However this must be done **all the way round** the rim. It is advisable to start at one point with both hands and then work, either all the way round the circumference in a circle, or if well used to the procedure, work both hands away from each other until they meet at the far side.



CLOSING WITH LIP TURNED UP

When sealing the containers with the lip turned up on the lid, fold the lip down and then repeat the closing process detailed above.

SPECIFICATION SHEET


SMALL CONTAINERS

No handles.


Supplied with WHITE lids.

RED, BLUE, GREEN & YELLOW
lids available in quantities of 100.

Please specify when ordering.

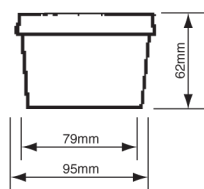


250ML
P40250

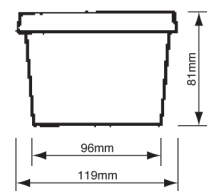


500ML
P40500

Net capacity underside of lid: 270ml
Gross capacity to rim: 300ml
Net capacity 10mm from top: 250ml


Net capacity underside of lid: 597ml
Gross capacity to rim: 639ml
Net capacity 10mm from top: 530ml

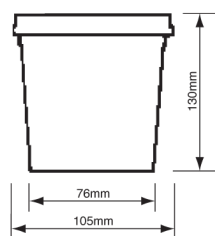

750ML
P40750

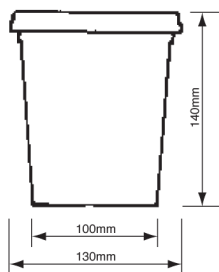

1000ML
P41000

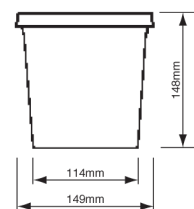

1.5 L
P41500

Net capacity underside of lid: 760ml
Gross capacity to rim: 830ml
Net capacity 10mm from top: 745ml

Net capacity underside of lid: 1170ml
Gross capacity to rim: 1196ml
Net capacity 10mm from top: 1122ml

Net capacity underside of lid: 1570ml
Gross capacity to rim: 1690ml
Net capacity 10mm from top: 1470ml


LARGE CONTAINERS

No handles.


Supplied with WHITE lids.

RED, BLUE, GREEN & YELLOW
lids available.


Please specify when ordering.

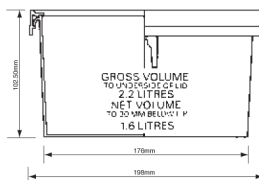


1.8 L
P41800

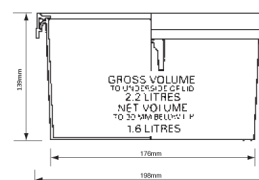


2.5 L
P42500

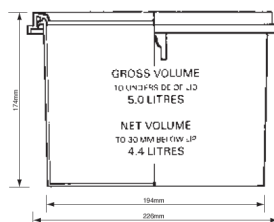


5.0 L
P45000

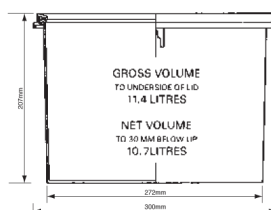


10.0 L
P410L

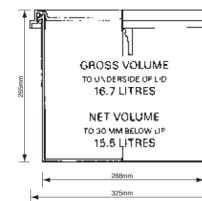

15.0 L
P415L

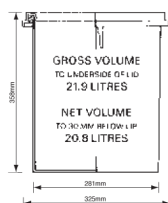

20.0 L
P420L

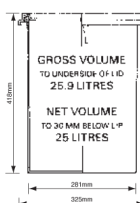

25.0 L
P425L

OTHER SIZES AVAILABLE

Please enquire about
other ranges of containers
in plastic and glass.

For further information or to request a sample for approval please contact our sales team.

SMART RELIABLE TRUSTED

WWW.SOLMEDIALTD.COM



TEL: +44 (0)844 8080 900

FAX: +44 (0) 844 8080 901

EMAIL: LABSUPPLIES@SOLMEDIALTD.COM